



ZETA POTENTIAL ANALYZER

## STABINO ZETA

**For the accurate determination of the zeta potential and colloidal stability, the STABINO ZETA is the first choice. It can replace the classical zeta potential measurement and is able to perform very fast titrations.**

Nowadays, particle surface charge and interface potentials, such as the zeta potential and the streaming potential, are widely used to characterize the stability of suspensions, emulsions and nanoparticles. These parameters have established as a typical measure representing the electrostatic repulsion between particles.

The STABINO ZETA has a high resolution and data point density, which allows very fast, precise and reproducible zeta potential measurements. It is possible to measure the zeta potential of particles in the range of 0.3 nm to 300  $\mu\text{m}$  in a concentration range of up to 40 volume percent. Regarding the optimized measurement technology, the STABINO ZETA can measure up to 5 parameters simultaneously and within a few seconds: Zeta potential, streaming potential, conductivity, pH value, and temperature. In combination with our unique NANOTRAC FLEX, the particle size can also be measured as a sixth parameter simultaneously in the same sample.




ZETA POTENTIAL ANALYZER STABINO ZETA

**FLAWLESS TITRATION**


The STABINO ZETA also has a built-in titration function in which all parameters are determined simultaneously at each titration dosing step. The determination of the isoelectric point is one of the titration options and is determined within a few minutes. Your titration options are:

**PH TITRATION**

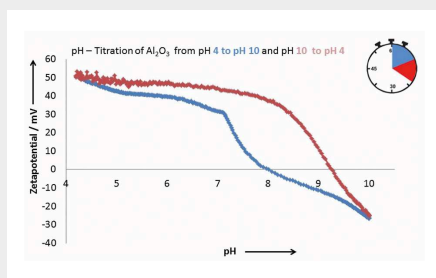


- | Determination of the isoelectric point
- | Stable pH ranges

**POLYELECTROLYTE TITRATION**

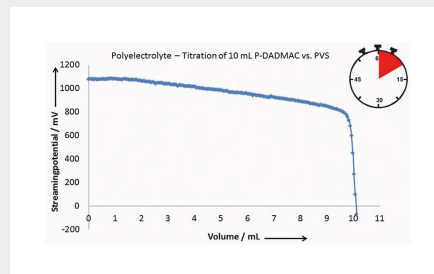

- | Statements about stability
- | Charge density
- | Dispersant optimization
- | Optimization in the formulation of your products

**TITRATION WITH SALTS**

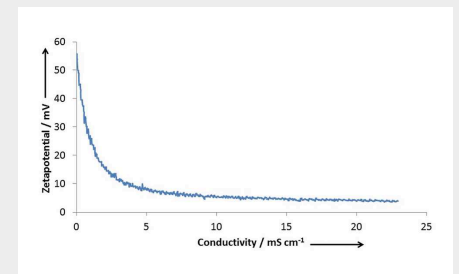

- | Zeta potential as a function of conductivity

ZETA POTENTIAL ANALYZER STABINO ZETA

## YOUR ADVANTAGES AT A GLANCE

### 5 MEASUREMENT PARAMETERS AT THE SAME TIME

To determine the quality of your samples, do you need more than just one measurement parameter? The STABINO ZETA provides you with information about the conductivity, zeta potential, streaming potential, temperature and pH of your sample with each measuring point.

### "MIX AND MEASURE" - AN ENORMOUS ADVANTAGE

Due to the continuous and rapid mixing of the sample and the titration solution, a charge titration is completed in minutes and additionally prevents sedimentation.

### MEASURING DURING TITRATION

With the STABINO ZETA software you can follow your entire titration or measurement in real time by means of the curve progression, because for each titrated drop you receive a measuring point with all 5 measuring parameters.

### ADJUSTED TITRATION SPEED

The titration speed of the STABINO ZETA can be adapted to the reaction speed of your sample. For this purpose, the software offers the possibility to define standard operating procedures (SOPs) as desired.

### FAST MEASUREMENT TIME

Most known analytical systems are based on the zeta potential of electrophoresis, where titrations are often too inaccurate and time-consuming. For high sample throughput and thus valuable time savings, the STABINO ZETA has been optimized so that parameters required for quality assurance, for example, can be determined within seconds. For a polyelectrolyte or pH titration, the STABINO ZETA requires only 5 - 15 minutes and can record several hundred data points.

### SIMPLE OPERATION

To focus only on the results, the software has been made as easy-to-use as possible. Just pour in 1 - 10 mL sample into the Teflon beaker of the STABINO ZETA, open the software and start the measurement.

ZETA POTENTIAL ANALYZER STABINO ZETA

## **NANOTRAC DUO – PRECISION IN PARTICLE AND ZETA POTENTIAL ANALYSIS**

The DUO unites two proven technologies – NANOTRAC FLEX and STABINO ZETA – into one complete characterization suite. This integration allows simultaneous measurement of size and zeta potential within the same sample, enhancing efficiency and providing comprehensive data.

Experience the DUO Advantage:

- | Comprehensive Analysis: Simultaneous size and stability measurements
- | Sample Preservation: No additional dilution required
- | Time Efficiency: Rapid, sequential measurements in one workflow
- | Data Correlation: Direct correlation between particle size distribution and zeta potential results
- | Flexible Use: Each instrument can be operated independently or seamlessly together as a single integrated solution



Applications:

- | Nanomaterials & Advanced Materials
- | Pharmaceutical Suspensions & Biotech Solutions
- | Coatings, Paints, and Pigments
- | Food, Beverages, and Nutraceuticals
- | Chemical Manufacturing & Polymers
- | Environmental Water Testing

ZETA POTENTIAL ANALYZER STABINO ZETA

**TOEBEHOREN**



Measuring cell 1 ml and  
3 ml with pestle



Measuring cell 10 ml -  
black -



Tempered measuring  
cell 0 - 90 °C



Piston set: 100  $\mu$  - 200  $\mu$   
- 400  $\mu$  - 1000  $\mu$  - 1200  $\mu$   
- 1500  $\mu$  - 2000  $\mu$  -  
conical

## ZETA POTENTIAL ANALYZER STABINO ZETA

### TYPISCHE TOEPASSINGEN

The STABINO ZETA is a highly versatile solution for rapid and reliable zeta potential and stability analyses. Designed to meet the demands of modern industries, it empowers users to optimize performance across a wide range of applications, including inks and pigments, ceramics, food and beverages, colloidal systems, polymers, microemulsions, cosmetics, battery slurries, chemicals, and carbon materials. Whether improving product quality, accelerating development, or ensuring process consistency, the STABINO ZETA delivers fast, actionable insights where they matter most.



*pigmenten / inkten*



*battery slurries*



*milieu*

- | pigmenten / inkten
- | keramiek
- | voedsel + dranken
- | colloidale suspensies
- | polymeren

- | micro-emulsies
- | cosmetica
- | battery slurries
- | chemicaliën
- | koolstof

- | milieu
- | zelfklevers
- | industriële mineralen
- | farmaceutica

... en meer!

Om de beste oplossing te vinden voor uw deeltjes-karakterisatie behoeften, kunt u onze toepassingsdatabase consulteren

ZETA POTENTIAL ANALYZER STABINO ZETA

**TECHNISCHE GEGEVENS**

<b>Methode</b>	Zeta streaming potential
<b>Berekeningsmodel</b>	none, as calibrated
<b>Metingshoek</b>	none, as mechanical measurement
<b>Size measurement</b>	yes (only in combination with NANOTRAC FLEX)
<b>Meetcel</b>	Teflon (10 ml, 3 ml, 1 ml)
<b>Zeta potentiaal analyse</b>	Ja
<b>Zeta streaming potential analysis</b>	Ja
<b>Zeta measurement range (charge)</b>	-3000 mV - +3000 mV
<b>Zeta meetbereik (grootte)</b>	0.3 nm - 300 µm
<b>Electroforetische mobiliteit</b>	Max. 14 (µm/s) / (V/cm)
<b>pH measurement</b>	Ja
<b>pH measurement range</b>	1 to 14
<b>Conductiviteitsmeting</b>	Ja
<b>Conductiviteitsbereik</b>	Up to 350 mS cm <sup>-1</sup>
<b>Temperatuursbereik</b>	0°C - 90°C
<b>Temperatuursnauwkeurigheid</b>	± 0.1°C
<b>Temperatuurscontrole</b>	Ja
<b>Titratie</b>	Ja
<b>Titration endpoints</b>	pH, zeta potential, conductivity, volume and time
<b>Reproduceerbaarheid (grootte)</b>	Refer to NANOTRAC FLEX
<b>Reproduceerbaarheid (zeta)</b>	2% with standard dispersion
<b>Monster volume zeta meting</b>	0.95 ml - 10 ml
<b>Monster concentratie</b>	Up to 40 %
<b>Meetvloeistoffen</b>	Water, polar organic solvents, acid and base
<b>Vochtigheid</b>	90 % niet-condenserend
<b>Afmetingen (B x H x D)</b>	180 x 300 x 260 mm

[www.microtrac.nl/stabino-zeta](http://www.microtrac.nl/stabino-zeta)